

Extraction is a process whereby the prepared shellfish tissue is converted into liquid containing released toxins.

Methanol Extraction Method



Equipment required to perform Methanol Extraction:

- Graduated cylinder 10ml*
- Transport tubes 5ml*
- Filters*
- 600 ml plastic beakers*
- Extraction solution bottle

Solutions required to perform Methanol Extraction:

- 100% Methanol
- Water

*Available to purchase from Jellett as individual components or as an extraction kit.

This extraction method can be used with the ASP Rapid tests.

Cleaning Up

Clean all equipment used in 5% household bleach and rinse well in tap water before re-using with another sample to avoid cross contamination.